

Calmodulin-Sepharose Beads

03/15

Store at 4°C. Do not freeze

Cat. No.: 7934-1	1ml
Cat. No.: 7934-5	5 ml
Cat. No.: 7934-10	10 ml

FORMULATION: Supplied as 50% aqueous slurry in 20 % Ethanol.

LIGAND DENSITY: ~ 1.5 mg Calmodulin protein per ml of drained resin/beads.

PREPARATION: Calmodulin-Sepharose Beads are prepared by covalently coupling recombinant human calmodulin (Cat # 7838) to 6% cross linked Sepharose beads.

INTRODUCTION: Calmodulin is a highly conserved regulatory eukaryotic protein involved in various cellular processes such as glycogen metabolism, cytoskeletal control, neurotransmission, phosphate activity and control of NAD/NADP. The diverse functions of calmodulin rely on its ability to interact with, and regulate the activities of large number of its target proteins/enzymes. Calmodulin binds proteins usually through its hydrophobic sites, which are exposed due to conformational changes in the presence of Ca^{2+} . However, some target proteins interact with Calmodulin in a calcium-independent manner. Calmodulin-Sepharose beads allow for specific enrichment of known or unknown calmodulin-binding proteins from biological samples. In addition, Calmodulin-Sepharose beads provide a convenient tool for affinity purification of calmodulin binding peptide (CBP) tagged recombinant proteins. Elution of CBP tagged proteins could be achieved by simply adding chelation ligand / removing Ca^{2+} .

APPLICATIONS:

- Single-step purification of native calmodulin-binding proteins.
- Tandem affinity purification (TAP) of protein complexes.
- Purification of CBP tagged proteins.
- Purification of calmodulin-regulated proteins from all eukaryotic cells.

RECOMMENDED BUFFERS:

Lysis / Binding / Wash buffer:

50 mM Tris Buffer, pH 7.4, 50-200 mM NaCl, 2 mM CaCl_2

- Avoid EDTA or EGTA in the buffers.
- Beads compatible with common reagents: 5 mM DTT, 10 mM β -mercaptoethanol, 0.1% TX-100/NP-40, and protease inhibitors.
- Increase the NaCl concentration for stringent washes if needed

Elution buffer:

Tris buffer pH 7.4, NaCl 50-200 mM, EGTA 5 mM

RECOMMENDED PROTOCOLS:

Batch Purification

1. Prepare samples in a Ca^{2+} containing buffer.
2. Wash beads with dH_2O , then equilibrate with binding buffer containing Ca^{2+} (For e.g., 50 mM Tris, pH 7.4, 150 mM NaCl, 2 mM CaCl_2).
3. Add the equilibrated beads directly to the sample lysate and incubate for several hours to overnight with gentle mixing (4°C).
4. Pour the beads-lysate mixture into a column, and wash with Ca^{2+} containing binding buffer, until no protein can be detected in the flow-through.
5. Elute with EGTA elution buffer (for e.g., 50 mM Tris, pH 7.4, 150 mM NaCl and 5 mM EGTA).
6. For Ca^{2+} -independent calmodulin binding proteins, glycine buffer (pH 2.0), urea elution buffer or boiling in SDS-loading buffer can be used.

Column (Gravity) purification

1. Prepare samples in a Ca^{2+} containing buffer.
2. Fill a column with the desired amount of Calmodulin-Sepharose beads and allow the beads to settle.
3. Wash the column with at least 5 column volumes of dH_2O .
4. Then wash the column with 5 column volumes of Ca^{2+} containing binding buffer.
5. Load the sample solution onto the washed column.
6. Collect the solution and repeat step 5 & 6 several times if necessary.
7. Wash the column with 5-10 column volumes of the binding buffer, until no protein can be detected in the flow-through, (increase the NaCl concentration if needed)
8. Add EDTA containing elution buffer to elute the bound proteins. Collect the eluents and store at appropriate temperatures.
9. For Ca^{2+} -independent calmodulin binding proteins, glycine buffer (pH 2.0) or urea elution buffer can be used.

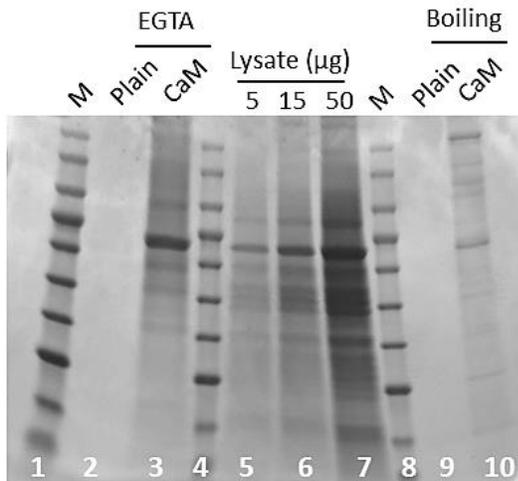


Figure: Enrichment of Calmodulin Binding Proteins (Calcium-dependent or -independent) using the Calmodulin-Sepharose: 293F cells were lysed with Lysis Buffer (50 mM Tris pH 7.4, 150 mM NaCl, 0.1% Triton X-100, 2 mM CaCl₂ and protease inhibitors). Clear lysate was collected after centrifugation at 16,000 g for 30 min at 4 °C. Calmodulin-Sepharose beads (CaM) and Plain beads (Plain) were washed with dH₂O, and equilibrated with Wash Buffer (50 mM Tris, pH 7.4, 150 mM NaCl, 2 mM CaCl₂). Approximately 25 mg clarified cell lysate (5.8 mg/ml) were incubated with 1 ml of Plain beads or CaM beads at 4 °C for 2 hours. After incubation, beads were extensively washed with Lysis Buffer, Wash Buffer, and finally with 500 mM NaCl in Wash Buffer. Beads were first eluted with 5 mM EGTA. Equal proportions (20%) of concentrated EDTA eluent were resolved on SDS-PAGE gel (EGTA, 2, 3). The proteins remaining bound to beads were eluted by boiling in 1x SDS loading buffer, and equal proportion of samples (5%) were used for SDS-PAGE (Boiling, 9, 10). Elution of proteins by EGTA and boiling is suggestive of Ca²⁺-dependent interaction (2, 3) and Ca²⁺-independent interaction (9, 10) respectively. This clearly indicates that large populations of proteins are captured by CaM beads (3 & 10), relative to the plain beads (2 & 9). Proteins eluted from CaM beads clearly display a distinct pattern from those in cell lysate (lane 5, 6, 7), supporting the enrichment of specific populations of proteins from lysates.

from those in cell lysate (lane 5, 6, 7), supporting the enrichment of specific populations of proteins from lysates.

RELATED PRODUCTS:

- Calmodulin, human recombinant (Cat No. 7838-500)
- Calmodulin, Bovine Brain (Cat. No. 7291-500, -1000)
- Active CAMK1b (Cat. No. 7729-5)
- Active CAMK1d (Cat. No. 7713-5)
- Active CAMK1G (Cat. No. 7736-5)
- Active CAMK4 (Cat. No. 7740-5)
- CaMKII Antibody (Cat. No. 3383-100)
- Phospho-CaMKII Antibody (Cat. No. 3384-100)
- Phospho-MARCKS Antibody (Cat. No. 3650-100)

FOR RESEARCH USE ONLY! Not to be used on humans.